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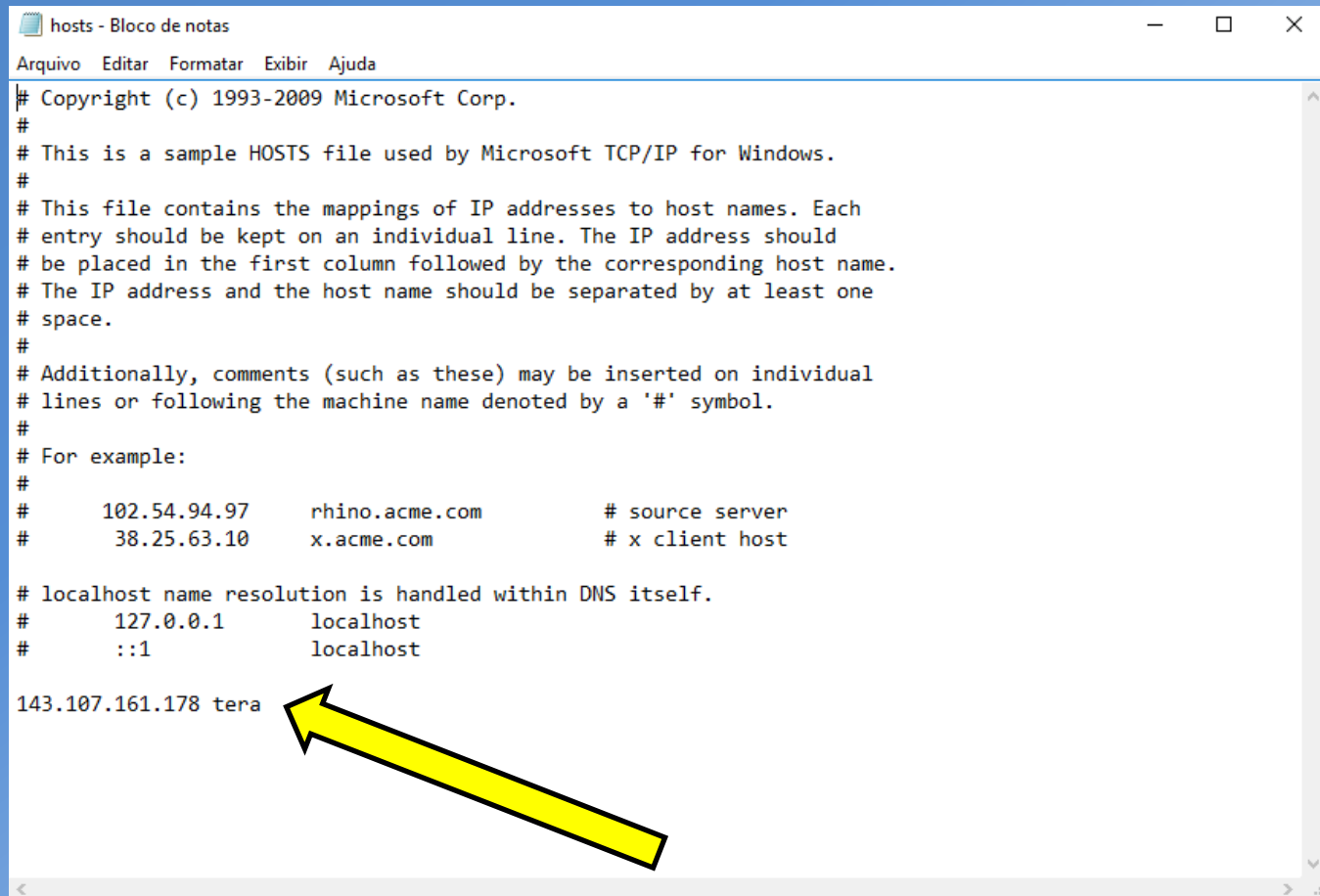
143.107.161.178 tera

```
hosts - Bloco de notas
Arquivo  Editar  Formatar  Exibir  Ajuda
# Copyright (c) 1993-2009 Microsoft Corp.
#
# This is a sample HOSTS file used by Microsoft TCP/IP for Windows.
#
# This file contains the mappings of IP addresses to host names. Each
# entry should be kept on an individual line. The IP address should
# be placed in the first column followed by the corresponding host name.
# The IP address and the host name should be separated by at least one
# space.
#
# Additionally, comments (such as these) may be inserted on individual
# lines or following the machine name denoted by a '#' symbol.
#
# For example:
#
#       102.54.94.97      rhino.acme.com          # source server
#       38.25.63.10     x.acme.com              # x client host
#
# localhost name resolution is handled within DNS itself.
#       127.0.0.1       localhost
#       ::1             localhost
143.107.161.178 tera
```

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
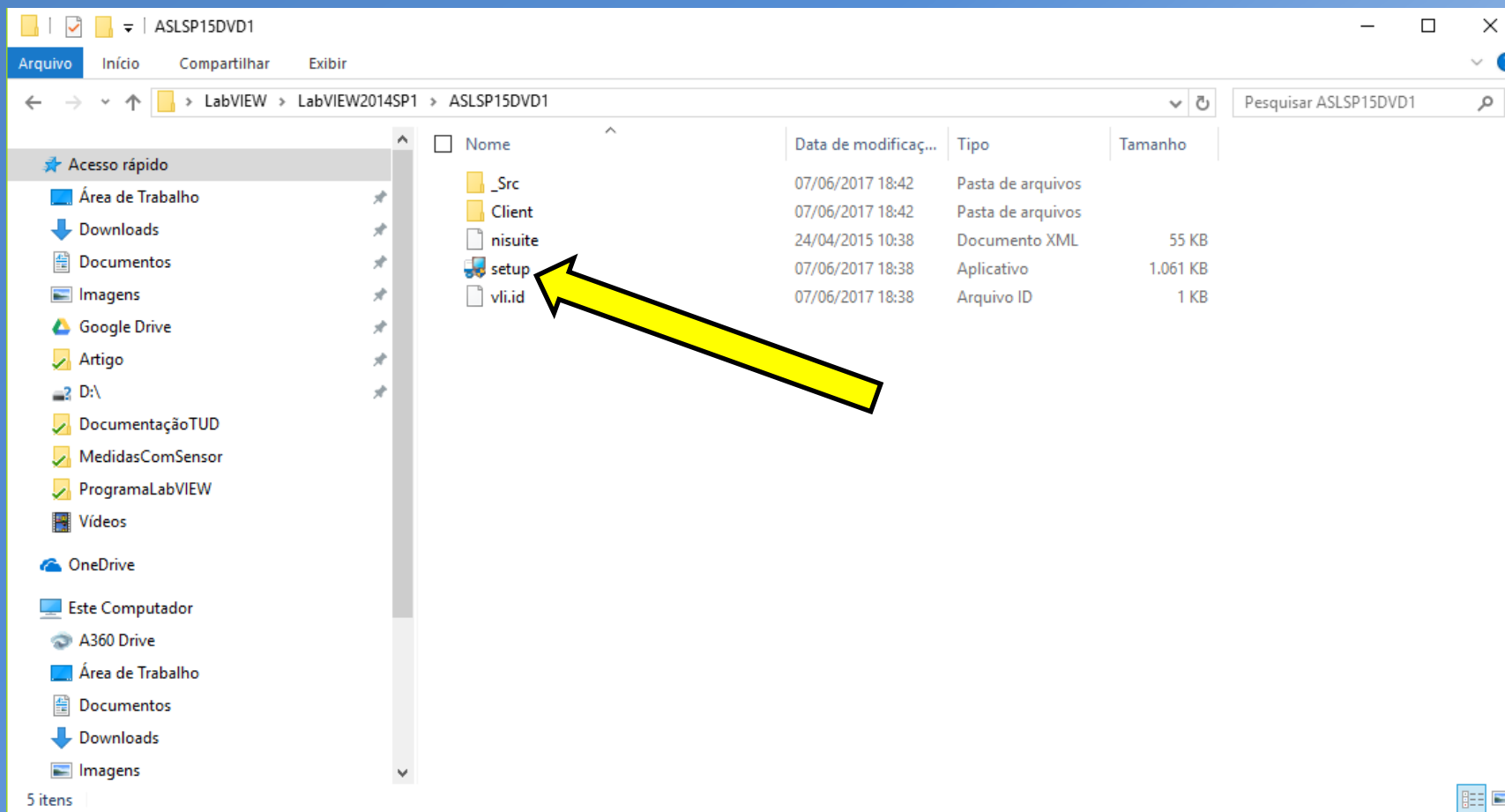
2) Caso necessite instalar os Devices Drivers, faça o download dos links:

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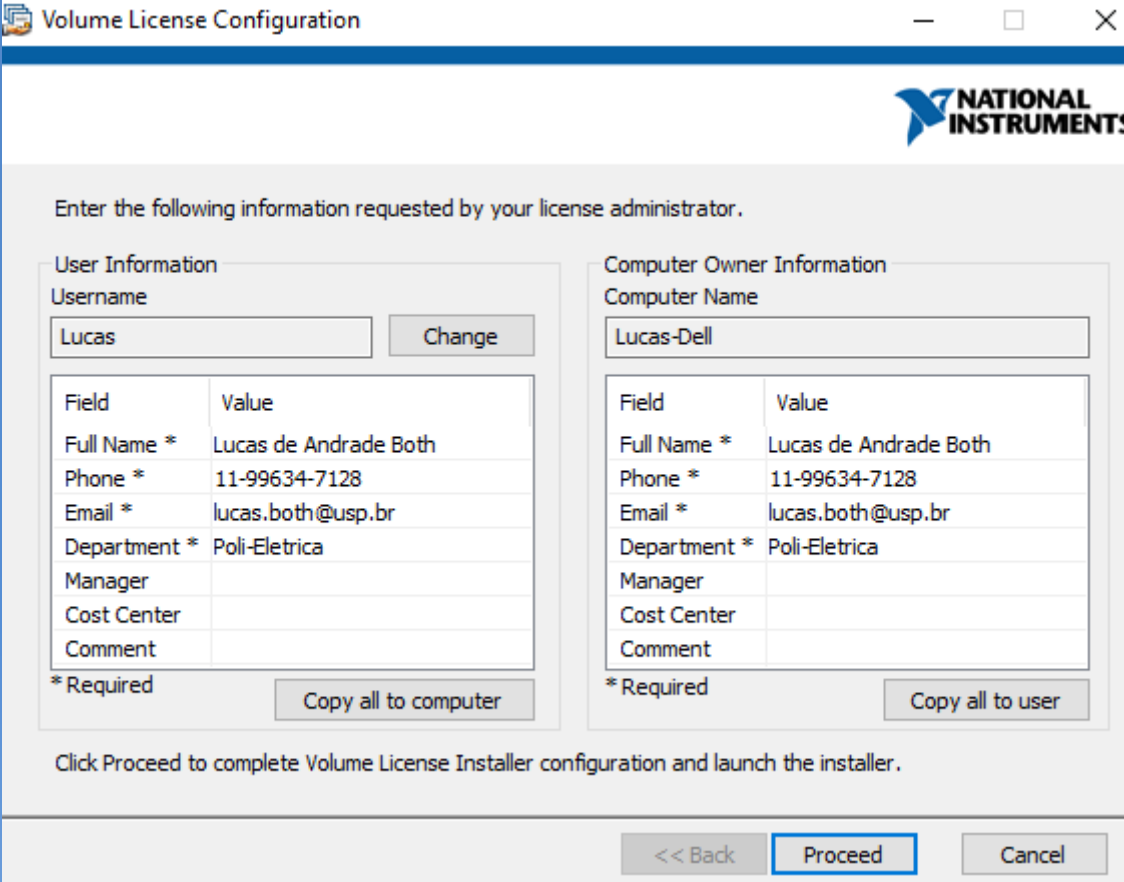The screenshot shows the 'Volume License Configuration' window with the 'Computer Owner Information' tab selected. The window contains two main sections: 'User Information' and 'Computer Owner Information'. Both sections have a 'Username' field and a table of user details. The 'Computer Owner Information' section also includes a 'Computer Name' field. At the bottom, there are navigation buttons: '<< Back', 'Proceed', and 'Cancel'. The 'Proceed' button is highlighted with a blue border.

Volume License Configuration

NATIONAL INSTRUMENTS

Enter the following information requested by your license administrator.

User Information

Username: Lucas

Field	Value
Full Name *	Lucas de Andrade Both
Phone *	11-99634-7128
Email *	lucas.both@usp.br
Department *	Poli-Eletrica
Manager	
Cost Center	
Comment	

*Required

Computer Owner Information

Computer Name: Lucas-Dell

Field	Value
Full Name *	Lucas de Andrade Both
Phone *	11-99634-7128
Email *	lucas.both@usp.br
Department *	Poli-Eletrica
Manager	
Cost Center	
Comment	

*Required

Click Proceed to complete Volume License Installer configuration and launch the installer.

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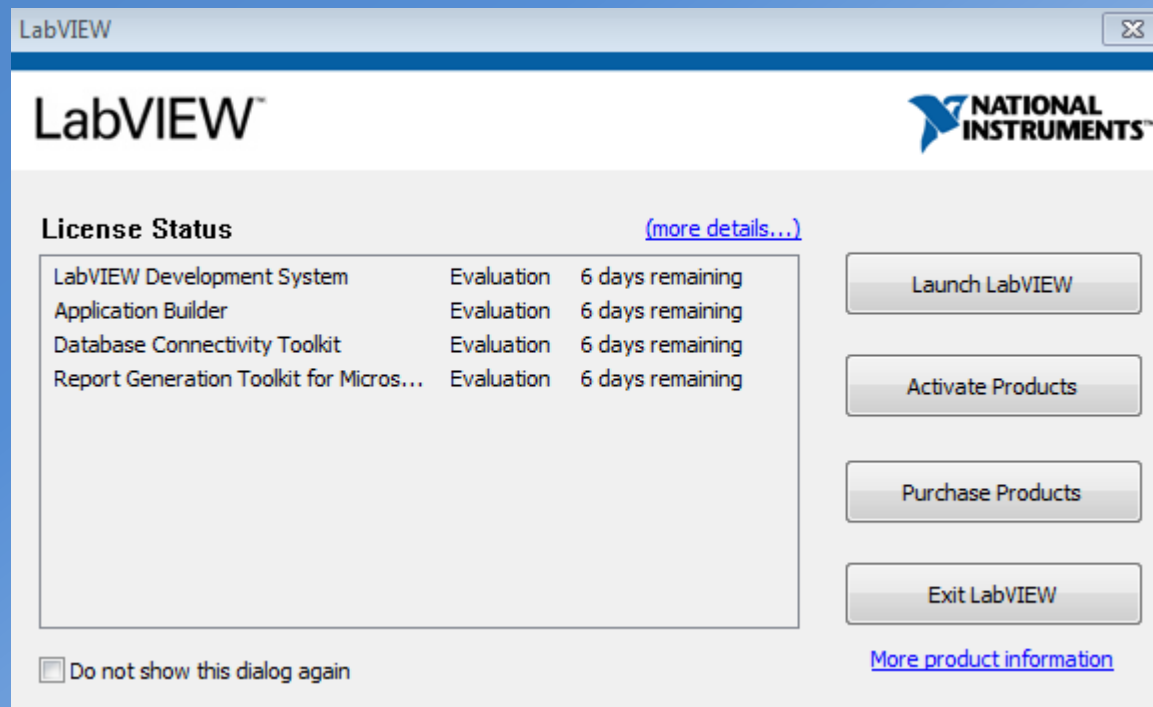
NI-VISA 15.0.1 - Pharlap ETS, NI Linux Real Time, VxWorks, Windows 10/8/8.1/7/Vista/XP (SP3) 32-bit/Server 2012 R2 64-bit/Server 2008 R2 64-bit/Server 2003 R2 32-bit

60 Ratings | 3.90 out of 5 |  Print

Available Downloads:

NI Download Manager (Recommended)

Non-Packaged Software Download: [NIVISA1501full_downloader.exe](#) (636.19 MB)
Checksum (MD5): 968498020cd9865291886c12a1b1344f

To get started:

- Click the **NI Downloader** link above.
- Your browser will download the **NI Downloader** launcher for your software (~150kb).
- Once **NI Downloader** launcher has been downloaded, launch the executable.
- The **NI Downloader** will automatically initiate the download of your software's standalone installer.
- When the download of the standalone installer is complete, launch it and follow the onscreen prompts to complete the installation of your software.

Note: Features of the NI Downloader include the ability to pause/resume download, or automatically resume download if unintentionally interrupted.

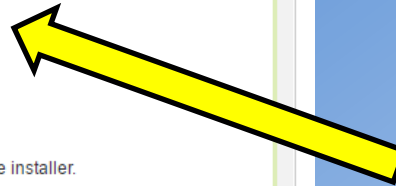
Browser Download

Browser Download: [NIVISA1501full.exe](#) (636.19 MB).
Checksum (MD5): 968498020cd9865291886c12a1b1344f

To get started:

- Click the **Standard Download** link above.
- Your browser will begin downloading the standalone installer for your software.
- Once the standalone installer has been downloaded, launch the executable and follow the onscreen prompts to complete the installation of your software.

Note: Browser managed downloads can be unintentionally interrupted due to dropped connectivity and do not provide the ability to "pause and resume".

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Drivers and Updates

Please fill out the information and click continue to retrieve your requested content.

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